

INDIRECT IMMUNOFLUORESCENCE

.... Fix cells.

Block 1 hour in 1:20 goat serum in 1X PBS.

Wash 5 mins with 1X PBS.

Add primary antibody in PBS in proper dilution.

Incubate 1 hour at RT.

Wash 3X with 1XPBS, once quickly, and twice for at least 20 minutes.

Add secondary antibody in PBS in proper dilution.

Incubate 1 hour at RT or overnight in the cold room.

Wash 3X with 1XPBS, once quickly, and twice for at least 20 minutes.

Mount slides with 50% glycerol + 10 mg/mL DABCO and store in the dark.